Item Check List 附件清单



Monitor 显示器




0.5 meter HDMI cable 0.5 米 HDMI 线



HDMI L shaped adapter HDMI 直角转接头



Y power splitter cable 一分二 DC 电源线

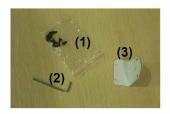

- (1) 12 Hexagon screws
- (2) L-shaped hex wrench
- (3) Brackets
 - (1) 螺丝12颗
 - (2) L型六角扳手
 - (3) 支架2个

Monitor Installation Instruction 显示器安装说明




- 1. Find the screws, L-shaped hex wrench and two white colored brackets that came with the monitor.
- 1. 找到附件中的螺丝和 L 型六角扳手和两个白色支架;



2. Place the monitor upside down on a clean table. As shown in the illustration, align two brackets to the lock pins on the back of the monitor.

Caution: Ensure there is <u>NO other object below the screen to avoid scratching</u> the screen surface when placing the monitor upside down on the table.

2. 显示屏平放在干净的桌子上,将支架按照如图方向放在显示屏支架孔。 注意: 当显示屏面朝下放置时,注意显示屏下面一定不要有其他物件,以免刮伤屏幕。

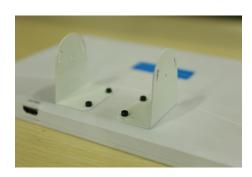

3. As shown in the illustration, use the screws to fix the brackets on the back of the monitor.

用螺丝将支架锁在显示器后面,四颗螺丝都锁好以后效果图如下:



4. Attach the TrueChrome camera to the brackets.

As shown in the illustration, face the label side toward you and place the camera between the brackets.

4. 安装 Truechrome II 相机: Truechrome II 相机有标签的一面在手心朝外,放入支架中;



- 5. Align the screw pins on the camera case to the lock pins on the brackets and use the screws to fix the camera to the brackets.
- Screw in a screw to the middle pin a bit first, but do not screw it in fully to allow to adjust the camera location angle.
- Adjust the camera position to make the screw pins on the camera case align to


the other two lock pins on the brackets and then tighten all the screws.

5. 螺丝孔对准后,先定位中间的螺丝(锁进去一点点,先不锁紧),再调整俯仰角度后再定位 另外两颗螺丝,三颗螺丝都定位好后再一起锁紧;



- 6. Screw in the screws on the other side of the bracket same as Step 5.
- 6. 另外一边的安装方法同第五步。

Cable Connection Instruction 接线说明

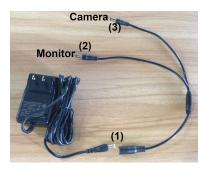

器,一路接相机。

- 1. Power Cable Connection
- As shown in the illustration, connect the 12V/2A AC/DC power adapter to the Y power splitter cable.
- Connect Y power splitter cable to the camera and the monitor.
- 1. 电源线连接方法

12V/2A 电源接一分二 DC 电源线,分为两路,一路接显示



2. HDMI Cable Connection

- Attach the HDMI L-shaped Adapter to the 0.5 meter HDMI cable.
- Connect the HDMI cable to the camera and Monitor
- 2. HDMI 线连接

将转接头连接在 0.5 米的 HDMI 线上。

HDMI 线一端接相机 HDMI 接口,一端接显示器 HDMI 接口。

